

SDS PAGE		
10% gels		
running	2 gels	1 gel
40 % acrylamide mix	5.000	2.500
1.5 M Tris HCl pH 8.7	5.000	2.500
SDW	9.600	4.800
10% SDS	0.200	0.100
10% APS	0.200	0.100
TEMED	0.020	0.010
<i>total</i>	<i>20.020</i>	<i>10.010</i>
stack		
40 % acrylamide mix	1.250	0.625
1M Tris HCl pH 6.8	1.250	0.625
SDW	7.000	3.500
10% SDS	0.100	0.050
0.1% BPB	0.300	0.150
10% APS	0.100	0.050
TEMED	0.050	0.025
	<i>10.050</i>	<i>5.025</i>
Pour away bubbles before running		
Run at 25 mA per gel		
Blot at 0.25 A for 60 min (two gels)		